## XEP 程式之掃描影像存檔基本設定

by 朱華偉 Sep,2005

 本實驗室 XE120 使用 XEP 軟體進行實驗掃描及操控功能。目前設定為一開 啟 XEP 即會自動跳出 Session Manager 視窗,提供使用者進行設定實驗資 料存檔路徑。所有 XE120 使用者,務必正確設定存檔路徑,以利 XE120 電 **腦管理者進行檔案管理。違規者實驗檔案可能會遭到刪除。**

Create Session	
Session Name	
Location	
0	
Upen previous session recently	used
)pen Session	
093005	
remanning	
romopping	
spmdata	
spmdata	
spmdata	
Location	
Location	

2. 欲設定新路徑請取消 "Open previous session recently used", 改用"…" 設定 欲儲存之資料夾路徑。設定完畢,按OK 按鈕。

Create Session	
Session Name	
093005	
Location	
C:\spmdata\test	
Open previous session rec Open Session 093005	ently used
Open previous session rec Open Session 093005 remapping spmdata	ently used
Open previous session rec Open Session 093005 remapping spmdata Location	ently used

3. XEP 採自動存檔檔名編排設定,使用者可利用工具列"Tools"下之 Preference 選項進行檔名編排設定。

Y XE Data Acquisition Program (PSIA) - [*XE-120* MFM XY:HIGH Z:HIGH CAN1								
🝸 File View Mode Setup	rools Help							
T 🗷 🖌 🔜 🗡	Layout Manager Session Manager	] 🔚 🥸						
📳 🐺 💾 🕰 🔇	Preferences	📲 💌 🔁 📕						
Scan Control	XEI (for Image Processing)							
□ Repeat ▼ Two way   X ○ X ○ Slow scan reverse	OY	Group ID : Not generated / Read 2µm /div						
Slope 0.000 Scan OFF	age -	-						

4 選取 Preference 視窗之 "Filename ",即可選定檔案自動儲存之檔名編排方式,按 "?" 即會列出所有編排方式的寫法及範例。請使用者自行參考使用。

n	Preferences			
Zoom 100 % V	Display Filename Image Motor Control Startup	Output filename ge Filename Format Filename Ex.	eneration rule [%1%N <date><sourcename><gennumb 050930Topography111</gennumb </sourcename></date>	er>
Output Filename           File name format striit           ## = <gennumber>           %1 = <date> (gymma           %2 = <time> (hhmm           %D = <date> (gymma           %D = <date> (gymma           %D = <date> (gymma           %D = <groupnumbe< td="">           %H = <hour> (D1, 12)           %H = <mont< td="">           %N = <mont< td="">           %N = <sourcename< td="">           %S = <sec> (D5, 30, %           %Y = <year> (D3, 04           Examples:           "PSIA_%1%4" creats           GenNumber means           Generation Number</year></sec></sourcename<></mont<></mont<></hour></groupnumbe<></date></date></date></time></date></gennumber>	Generation Help g contains the following identifie (001, 020, 055, 100,) id/030430, 031230,) ss/130401, 020530,) onth> >> (001, 002, 003,) ,18,) y5,] y4, 12,} > (Topography, Z Detector,) 55,] es the output filename "/SIA_ <c he output filename "/SIA_<c he output filename "/SIA_<c< td=""><th>ars Date&gt;%4<gennumber>.tiff". nNumber%5.tiff". inNumber%5.tiff". i the output file name.</gennumber></th><td>#", Err. Sig. Z Scanner</td><th>Cancel Apply 0000 X=+00.0000 Y=+00.000 X Z Stage Lift Z</th></c<></c </c </c </c </c </c 	ars Date>%4 <gennumber>.tiff". nNumber%5.tiff". inNumber%5.tiff". i the output file name.</gennumber>	#", Err. Sig. Z Scanner	Cancel Apply 0000 X=+00.0000 Y=+00.000 X Z Stage Lift Z